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- GB** Detach a swab from block (original is fully packed by aluminium packaging) and remove one side of the packaging by holding on to the other side.  
Note: Do not throw out this packing. It will be used later.

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- GB** Swab required test site using the open side.

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- GB** In case of positive (existence of protein residues greater than detection limit of 20µg), the swab will change to green color within seconds for high protein concentration and take up to 2 minutes for low protein concentrations.

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- GB** Another test site can now be tested with the unused swab on the other side by placing the cover back on the already used swab.

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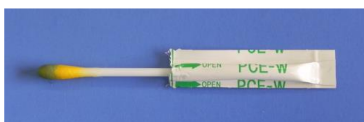
18 months after manufacturing

**Precautions**

- After rubbing check point with Pro-Check, its pigment may remain at the surface. Wipe out the pigment by washing with water or cleaning with alcohol. The pigment is harmless.
- If the test point is asperity, the pigment may remain at concave side. In such a case, cleaning can be carried out by Chlorine detergent or acetic acid.
- A part of alcohol might be evaporated from this product during storage. In this case, swab will change to red from yellow. You will be able to use the swab after moistening it with alcohol for disinfection.
- The swabs must be used within an hour of opening.

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